Référence **215670**

## **Kit Vérification MD100 MD200**

### **Verification Standard kit Photometer MD100 MD200**

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*The verification standard kit for the MD 100 / MD 200 is designed to assure the user of the accuracy and the reliability of the results related to the integrated wave lengths.*

*The kit contains one zero standard, 6 different vials for checking 6 different wave lengths and allows checking the complete range of MD 100 photometers.*

*The shelf life of the verification standard kit is two years from the date of production, provided that storage and use are in accordance with the instructions provided.*

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## Absorbance measurement using a Photometer MD 100 and MD 200

To measure Absorbance, press the [Zero/Test] key and hold. Turn the instrument on by pressing the [On/Off] key. Once three decimal points appear in the display, release the [Zero/Test] key.

The display temporarily shows: **AbS** followed by the wavelength to be measured (e.g. **A530** for 530 nm).

For instruments with more than one LED, the desired wavelength can be selected by scrolling through the list using the [Mode] key.

Select the vial marked with „Zero“, place it in the sample chamber and cover the vial with the light shield. Press the [Zero/Test] key and allow instrument to zero. Remove the 'Zero' vial.

Next, select the vial containing the reference standard marked with the appropriate wavelength (e.g. A530), swirl gently several times, and place it in the sample chamber. Cover the vial with the light shield provided and press the [Zero/Test] key.

Allow unit to take measurement and compare the reading on the display with the values specified in the reference standard certificate, taking the tolerances into account (see also certification notes & directions).